

**GENERAL DESCRIPTION**

The Septihol™ WFI Sterile Alcohol Wipes combine pre-saturated knitted polyester wipes with a ready-to-use (RTU) solution of 70% (v/v) USP Isopropanol and 30% (v/v) USP Water for Injection (WFI). This right-sized packaging system provides a convenient method to introduce low particulate wipes to your production environment while reducing product waste. Each package is double bagged, and gamma irradiated for use in aseptic areas. The outer pouch has a resealable label/closure. Septihol WFI Sterile Alcohol Wipes are part of the Septihol family of products. Each product is designed to simplify your validation requirements while providing multiple packaging options.

The Septihol WFI Sterile Alcohol Wipes are intended for the cleaning and removal of residues from hard non-porous surfaces in cleanrooms or other critical environments such as floors, walls, counters, tools, carts, and other equipment.

**FEATURES**

**BENEFITS**

70% (v/v) USP Isopropanol and 30% USP WFI alcohol solution

Formulation easily evaporates

Right-sized packaging

Knitted polyester wipe with sealed edges

Double bagged and gamma irradiated, with every lot USP sterility tested

Manufactured under FDA cGMP, ISO 9001 and ISO 13485 compliant quality system

Industry best practice for cleaning and rinsing

Ideal for use on critical surfaces

Reduced waste of wipes

Low particulate wipes to achieve compliance with current Annex 1 contamination control strategies

Facilitates introduction into critical areas

Supports quality assurance and quality control

**PHYSICAL PROPERTIES**

**Septihol WFI Sterile Alcohol Wipes**

Size ..... 12x12 in (30.5 x 30.5 cm), 15 sheets/pack

Color ..... White

Material ..... 100% knitted polyester, sealed edge

Fibers ( $\geq 100\mu\text{m}/\text{cm}^2$ ) .....  $\leq 0.20$ , unsaturated

Particles ( $\geq 0.5\mu\text{m}$ ) ..... (Biaxial Shake x  $103/\text{cm}^2$ )  $\leq 1.0$ , unsaturated

Specific Gravity (25 °C [77 °F]) ..... 0.872 typical

Fill Volume (mL) ..... 427, typical

Sterility Test ..... Sterile, Current USP

**DIRECTIONS FOR USE**

The Septihol WFI Sterile Alcohol Wipes are designed to be used on hard non-porous surfaces in pharmaceutical, biotechnology, cosmetic, nutraceutical, medical device, university and government research, and laboratory animal research facilities that employ cleanrooms and other controlled environments.

Remove the double bag system. Ensure package opening is facing up. Using the resealable cover, remove wipes. Reseal once desired number of wipes have been removed. Using the wipes, wet surfaces thoroughly by spreading manually with appropriate application technique. Allow surfaces to air dry.

The Septihol WFI Sterile Alcohol Wipes are a complete product. Do not add other chemicals.

## STORAGE

---

Do not contaminate water, food, or feed by storage or disposal. Open dumping is prohibited. Do not reuse empty pouches or used wipes. This product should be stored in an area where it will not be exposed to extreme temperatures.

## DISPOSAL

---

Dispose wipes in an appropriate waste container. Dispose of in accordance with local, state, and federal regulations.

## SERVICE

---

### Sales

Service is one of the most important ways to verify consistent quality of the facility's performance and operation. A tailored service program by STERIS provides effective, trouble-free operations.

### Technical

STERIS is pleased to provide a completely staffed and equipped technical service laboratory capable of performing needed tests and providing both telephone and on-site assistance when needed. More details on how this service can benefit a facility's particular situation can be provided upon request.

## PRECAUTIONS

---

Information concerning human and environmental exposure may be reviewed on the Safety Data Sheet (SDS) for the product. For additional information regarding incidents involving human and environmental exposure, call 1-314-535-1395.

---

### For further information, please contact:



STERIS Corporation  
5960 Heisley Road  
Mentor, OH 44060-1834 • USA  
440-354-2600 • 800-444-9009  
[www.STERISLifeSciences.com](http://www.STERISLifeSciences.com)